
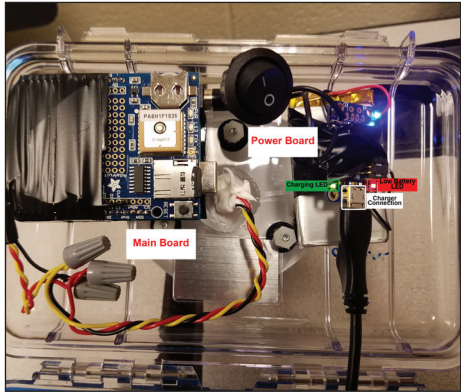


## Sensor Operation and Charging

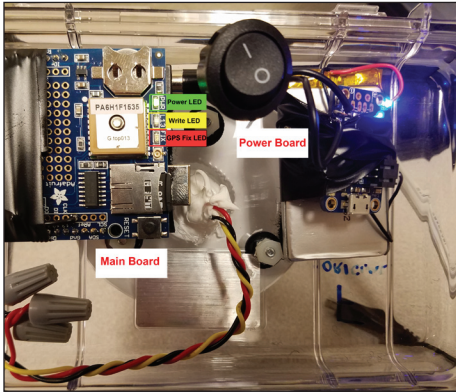

## Sensor Operation and Charging

### To Charge:

- Flip power switch to “off” position.
- Connect charger to power unit.
- Make sure green charging LED is lit.
- When charging is complete, unplug the charging cable.
- When battery is low, the red low battery LED will light up.

### To Operate:

- Disconnect charger and flip power switch to “on” position.
- Make sure the green power LED is lit.
- Make sure that the lip of the lid is clear of obstruction; close and latch lid.
- The yellow write cycle LED flashes whenever data are being written.